

# Passive Diffuser Cell

Model: 7011



## Application

This passive diffusion cell for tritium is used to determine the total tritium in air based upon taking a continuous air sample where the vial is located.

## Features

- Simple
- Inexpensive
- Effective
- Accurate
- Reliable

## Description

The Passive Diffusion Cell is a rudimentary but effective apparatus that should be used in all tritium laboratories to measure airborne tritium activity. It comprises a vial filled with water and fitted with 5 micron diffuser plates installed inside the 1/2" stainless steel inlet tube inserted through the lid.

The vial is held in position in a yellow stand, to which radiation sticker clearly identifying it as a tritium instrument is attached. The stand is intended for a table or shelf, or can be mounted on a wall.

As the Passive Diffuser Cell sits in the atmosphere of the laboratory, it absorbs tritium from the air, and the uptake of tritium has a relationship to the amount of tritium present in the room atmosphere. The cell contents are measured weekly in a scintillation counter. With several of these inexpensive passive diffusion cells placed selectively around a laboratory, convenient mapping of the total tritium presence in the atmosphere is recorded. Readings obtained confirm values from other often more sophisticated but less reliable instruments, confirming total tritium presence and potential exposure by personnel. After use, the vial is cleaned and refilled with water in preparation for the next measurement.

# Specifications

Vial Material	Plastic
Vial Size	5 cc
Holder Size	5" x 5" x 6" high
Holder Material	Powder painted carbon steel

Agent Stamp / Business Card